



NOPA

NOPA

The NOPA bag is a sterile solution specially designed for sampling in the oilseed sector. It ensures reliable results while meeting the rigorous standards set by the National Oilseed Processors Association (NOPA). For decades, NOPA has established stringent standards and advocated for the oilseed processing industry's interests. Featuring a writing area, the NOPA bag guarantees optimal traceability.

It is the indispensable tool for the oilseed processing industry!

Features and Benefits

- Made from FDA-compliant material
- Meets NOPA and COPA standards
- Suitable for oilseeds
- Leak-proof and airtight
- Can be used in the homogenizer
- Guaranteed sterility
- Free from RNase, DNase, and pyrogens




Inspired by your challenges

LABPLAS

VISIT LABPLAS.COM FOR ADDITIONAL INFORMATION ABOUT OUR PRODUCTS | Tel: 1.450.649.7343



NOPA BAG INSTRUCTION

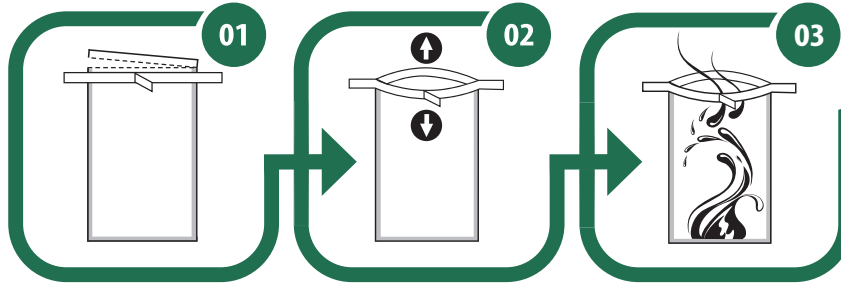

01
Tear off the top of the bag
along the perforations

02
Use pull-tabs
to open the bag

03
Aseptically insert the
sample in the bag



04
Pull the wire-ends on
each side of the bag

05
Twirl the bag around the
wires, at least 3 times,
to close the bag

06
Fold both wire-ends onto
the bag to provide airtight
and leak-proof closure



Inspired by your challenges

LABPLAS

VISIT LABPLAS.COM FOR ADDITIONAL INFORMATION ABOUT OUR PRODUCTS | Tel: 1.450.649.7343

